Quick Start

FTDI CDM Drivers

Extract the FTDI CDM Drivers.

FTDI CDM Drivers

Extract Cancel

To continue, accept the following license agreement. To read the entitiagreement, use the scroll bar or press the Page Down key.

IMPORTANT NOTICE: PLEASE READ CAREFULLY BEFORE INSTALLING THE RELEVANT SOFTWARE:

This licence agreement (Licence) is a legal agreement between you (Licensee or you) and future Technology Devices International Lin of 2 Seaward Place, Centurion Business Park, Glasgow (641 HH, Scotland (UK Company Number SC136540) (Licensor or we) for us

install

99 %

Exit

< Back Next > Cancel

You are only a few steps away from installing and hence, using a PhysLogger. Download the Setup.

Additional Installations

While installing,

PhysLogger Desktop

App requires some

download with it.


additional files to be



Oosain Package Man

Downloads

Features

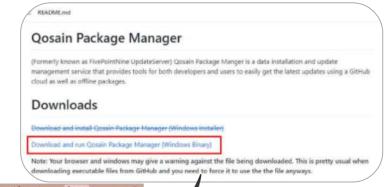

Windows protected your PC

Your browser may warn you against downloading this executable, but you need to ignore this message.

Run anyway Don't run

Installation Location

When warned by MS Defender SmartScreen. select More information and then click on Keep anyway.

Initiate Installation.



package manage

Once

For the Device

Driver Installation Wizard, select "I accept the agreement" and click Next.

PhysLogger Desktop

Overall

About the project The first stable release of PhysLogger Deskton

Back Next

Specify the directory where the

Recommended is the default one.

PhysLogger Desktop App

package will be installed.

Last Updated: 2022-Mar-17 05:39:54

PhysLogger I

The package would require 198.73 MBs on your computer

Last Lindated: 2022-Mar-17 05:39:54

out the project The first stable release of PhysLogger Desktop

Allow all required files and updates to be installed. Wait till the installation is 100% complete.

install PhysLogger Desktop Publisher: Qosain Scientific Last Updated: 2022-Mar-17 05:39:54 out the project. The first stable release of Physl on Click 'Extract' to unpack version 2.12.28.2 of FTDI's Windows driver package and launch the installer. FT232 drivers are required to run PhysLogger. Even if your windows automatically Introsoft DirectX (Web Installtion)
For the best viewing experience, PhysLogger requires the latest version of DirectX I Irdu Naetalaan Fonte Back Next Exit nstalling Microsoft(R) DirectX(R

DirectX

During the installation of MS DirectX when asked, uncheck the installation of Bing Bar.

For the auxiliary installation of Pak Urdu Installer, skip restart now and finish installation first.

Up-to-the-minute headlines at the click of a buttor

Assessment and Physics Strategies to the Microsoft Service

You'll get future updates to Bing Bar and other Microsoft products from Microsoft Update. This software may also download and install some updates automatically.

< Back Next > Cancel



Allow all the necessary drivers to be installed.

Desktop Shortcut

Once installation has completed successfully, you will see this PhysLogger Desktop dialogue box and a desktop shortcut icon on your screen. It will also appear in the Start menu.



Try the Bing Bartoday

K Back Finish Cancel